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Determination of naphtha in oxygen-containing organic compounds by gas chromatography

Naphtha (Naphtha) also known crude gasoline: containing full and naphthalene alkane 55.4%, single ring alkane 30.3%, double 2.4% of alkanes, alkylbenzene 11.7%, 0.1% benzene, indene with 0.1%, the main raw material for ethylene production. In naphtha methyl tert butyl ether (MTBE), containing oxide belongs to the impurities, once cracking feedstock in the containing oxide content is higher than 1000 mg kg⁻¹, will lead to cracking of propylene and ethylene alcohol.

The method provides for the methyl tert butyl ether (MTBE), ethyl tert butyl ether (ETBE), tert amyl methyl ether (TAME), methanol, ethanol, organic containing oxygen compounds by gas chromatography method for the determination of naphtha.

This method is suitable for determination of naphtha of the oxygen-containing organic compounds determination range: 5mg/kg ~ 1000mg/kg.

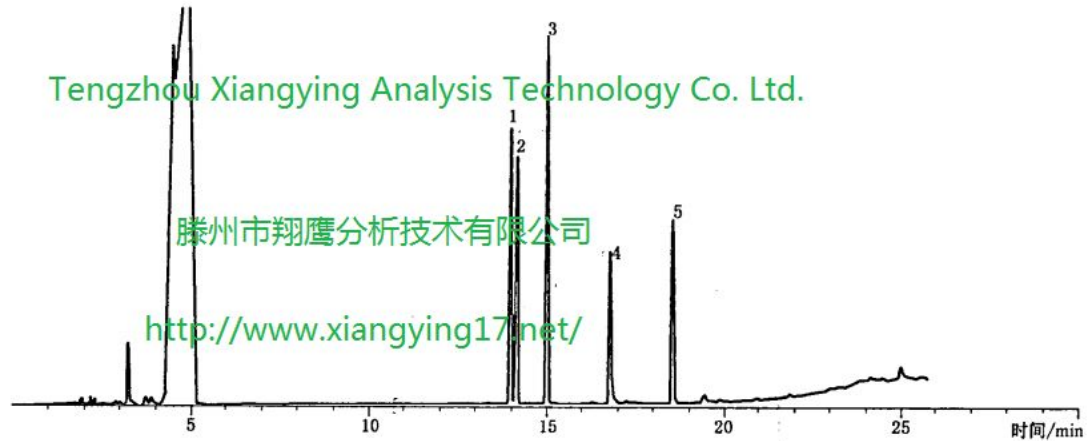
Tengzhou City Xiangying Analysis Technology Co., Ltd. the XiangyingGC7990plus gas chromatography determination in naphtha trace containing oxides, using a non tandem and independent separation of polar column and a polar column, completed within 30 min in naphtha trace containing oxides analysis. The system peak area reproducibility is good, and the RSD was less than 1.5%, quantitation limit and low limit of detection. The method is accurate and reliable.

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XiangyingGC7990plus features:

- 1 excellent performance, competitive price, to meet the customers' choice
 - 2 multi industry adaptability
 - 3 outstanding performance in all kinds of harsh environment
- INJ

A variety of sampling ports can be equipped with: packed column inlet, shunt / non shunt capillary inlet,

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