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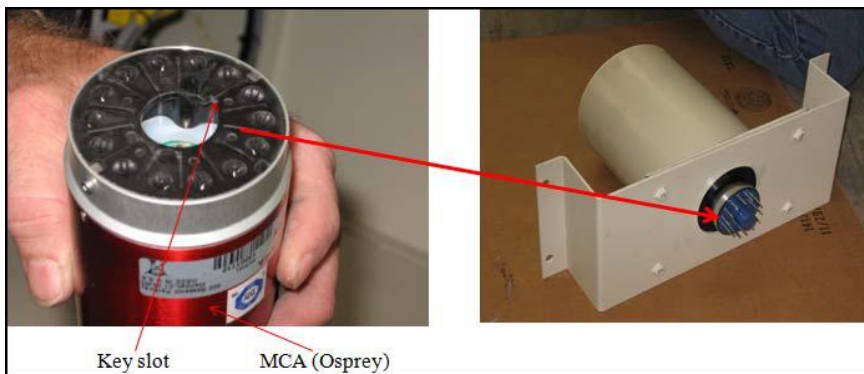
TEMCO, Inc.

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Instructions | Amplifier Settings | Calibration | Measurements

Belt Speed Guidelines

- 1) For total gamma measurements, set to a maximum of 1 feet/minute for open core; If the core is in inner barrel, set to a maximum of 0.5 feet/minute.
- 2) For spectral measurements, set the speed at a maximum of 0.5 ft/min.

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TEMCO, Inc.

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Instructions | **Amplifier Settings** | Calibration | Measurements

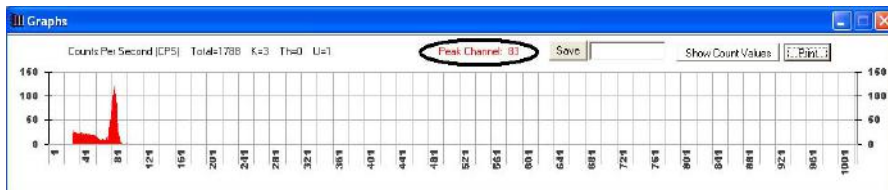
Acquisition Time (seconds)

Amplifier Coarse Gain

Amplifier Fine Gain

Amplifier High Voltage +

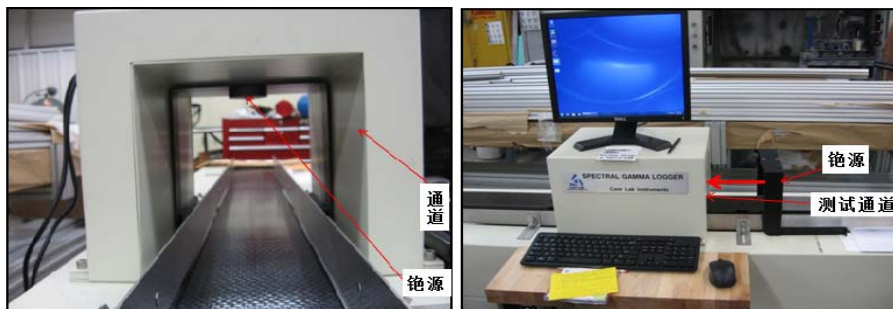
TEMCO, Inc.

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TEMCO, Inc.

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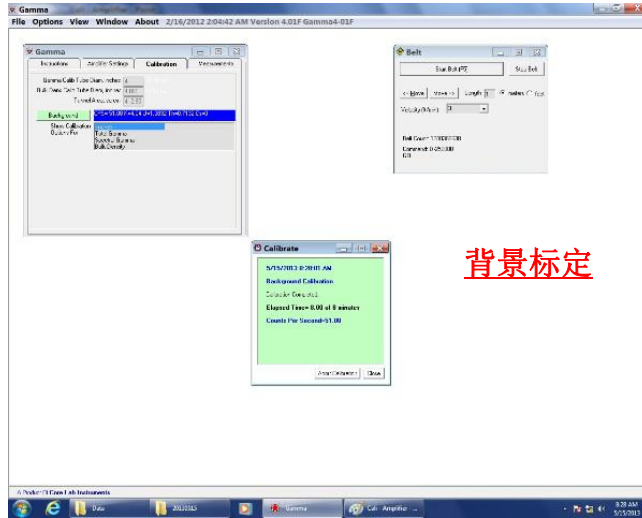
Instructions	Amplifier Settings	Calibration	Measurements
Calibration Tube Diameter <input type="text"/> <input type="text"/>			
Tunnel Area (sq cm) <input type="text"/>			
Background		CPS =	
Show Calibration Options For <input type="text"/>			
API Calibration		API Tube API <input type="text"/>	CPS =
Potassium Calibration		Potassium % <input type="text"/>	CPS =
Uranium Calibration		Uranium ppm <input type="text"/>	CPS =
Thorium Calibration		Thorium ppm <input type="text"/>	CPS =
Bulk Density Aluminium		Density, gm/cc <input type="text"/>	CPS =
Bulk Density Polycarb.		Density, gm/cc <input type="text"/>	CPS =

TEMCO, Inc.

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TEMCO, Inc.

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TEMCO, Inc.

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Gamma							
Instructions		Amplifier Settings		Calibration		Measurements	
Gamma Calib Tube Diam, inches	4						
Bulk Dens Calib Tube Diam, inches	4.000						
Tunnel Area, sq cm	412.90						
Background	CPS= 51.08 K=4.04 U=1.3832 Th=0.7132 Cs=0						
Show Calibration Options For	<none> Total Gamma Spectral Gamma Bulk Density					伽玛谱标定	
API Calib	API Tube API	200.00	CPS= 47.89				
Potassium (K) Calib	Potassium %	4.50	CPS= 1.8117%				
Uranium Calib	Uranium ppm	100.00	CPS= 3.9545				
Thorium Calib	Thorium ppm	100.00	CPS= 1.4979				

TEMCO, Inc.

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Gamma							
Instructions		Amplifier Settings		Calibration		Measurements	
Gamma Calib Tube Diam, inches	4						
Bulk Dens Calib Tube Diam, inches	4.000						
Tunnel Area, sq cm	412.90						
Background	CPS= 51.08 K=4.04 U=1.3832 Th=0.7132 Cs=0						
Show Calibration Options For	<none> Total Gamma Spectral Gamma Bulk Density					总密度标定	
Bulk Dens Aluminium	Density, gm/cc	2.700	CPS= 397				
Bulk Dens Polycarb.	Density, gm/cc	1.200	CPS= 978				

TEMCO, Inc.

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TEMCO, Inc.

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The screenshot shows the 'Gamma' software interface with the following data:

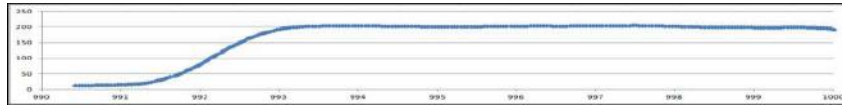
Field	Value
Job Name	Test - API200
Company	CNOOC SZ
Location	SZ
Well	NA
Operator	Henry
Bottom Depth	1000 Feet
Top Depth	998.5
Core Diameter	4 Inches
Cross Section	Unslabbed
Run Type	Total
Velocity (ft/min)	0.25
Depth At Detector	998.66
API	194

Additional settings visible: Slice Length (Inch)=1, Log Debug Information checked, File Name: Test - API200.

TEMCO, Inc.

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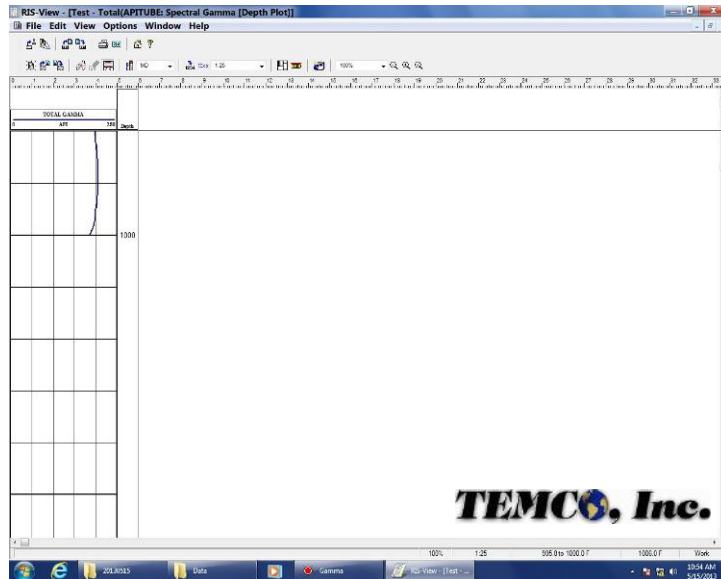
TEMCO, Inc.

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TEMCO, Inc.

**RADIOACTIVE MATERIAL
LIMITED QUANTITY - UN2910**

This package conforms to the conditions and limitations specified in 49 CFR 173.421; IATA Section 10.5.9.4; and Postal Publication No. 6 dated September 1983 for radioactive material, excepted package, limited quantity of material - UN2910.

The radiation exposure rate at any point on the external surface of this package does not exceed 0.5 millirem per hour.

No other labels are required - or authorized.

The radioactive material contained in this package is exempt from USNRC and Agreement State licensing requirements.

"Radioactive Material - Not for human use- introduction into foods, beverages, cosmetics, drugs, or medicinals or products manufactured for commercial distribution is prohibited - exempt quantities should not be combined" (from USNRC regulation 10 CFR 32.19d).

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Cochi (Beijing) Technology Limited